


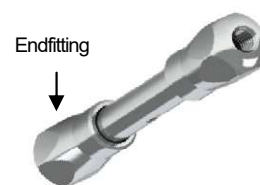


【Instructions For Column Connection YMC-Triart Metal Free Column】

To ensure optimal performance of the column, please be careful with the column connection.

- The column endfitting (see picture in right) is Parker style connection. Connect the column by hand-tightening. Using a wrench for column connection may damage the column.
- Make sure to hold the endfitting when connecting or disconnecting the column. Leakage may occur if the endfitting is loosened.
- A ferrule-integrated type connector may damage a frit in the column when it is over-tightened. In order to avoid such trouble, use of a connector which does not have a ferrule or a ferrule replaceable connector is recommended (See below for our recommendation).



Recommendation	◎	○	
Ferrule	NO	REPLACEABLE	
Product	MarvelX	Handy connector 2	Hand-tight EXP [®] fitting
Manufacturer	IDEX Health & Science LLC	YMC	Optimize Technologies, Inc.
			

*MarvelX is a registered trademark of IDEX Health & Science LLC

*EXP is a registered trademark of Optimize Technologies, Inc.

Please contact your local YMC office for more details.

YMC CO., LTD.

YMC Karasuma-Gojo Bldg., 284 Daigo-cho, Karasuma Nishiiru Gojo-dori, Shimogyo-ku, Kyoto 600-8106, Japan

TEL +81-75-342-4515 FAX +81-75-342-4550

URL:www.ymc.co.jp E-Mail:sales@ymc.co.jp

(UM170410AE)